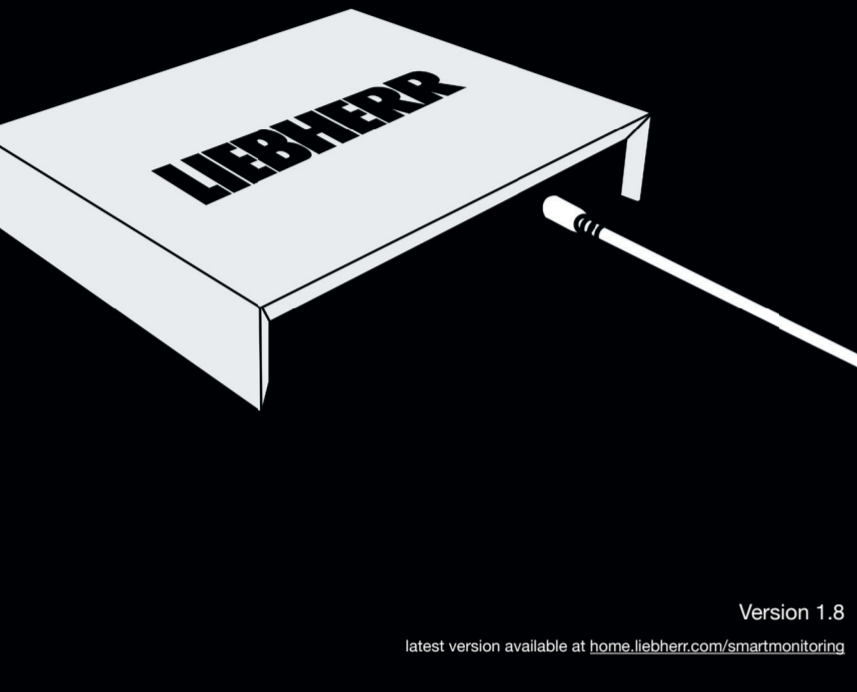
# SmartCoolingHub

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Version 1.8

latest version available at [home.liebherr.com/smartmonitoring](http://home.liebherr.com/smartmonitoring)

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For more information and other accessories visit the SmartMonitoring website: [home.liebherr.com/smartmonitoring](http://home.liebherr.com/smartmonitoring) or contact your dealer.

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<b>EN</b>	<b>Start the set-up at the following address:</b> <a href="http://smartmonitoring.liebherr.com/setup">smartmonitoring.liebherr.com/setup</a> or scan this QR-Code:

You will also find the following there:

- Set-up videos
- Appliance instruction manual
- SmartMonitoring instruction manual

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<b>EN</b>	Off	SmartCoolingHub is switched off
	Flashing green and red	SmartCoolingHub is starting up
	Flashing green rapidly	Ready for set-up
	Flashing green slowly	SmartCoolingHub is being set up
	Constant green	System is operating normally
	Constant red	Independent reboot of the SmartCoolingHub (may take several minutes).
	Flashing red rapidly	Network error (no IP address received or no Internet connection available)

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For more information, refer to the SmartMonitoring instruction manual.

- DHCP mode active (automatic IP address assignment)
- The MAC address of the SmartCoolingHub must be shared in the network (see label on the back of the SmartCoolingHub if a MAC filter is used in your network).
- An Internet-enabled device with a current browser version and a minimum resolution of 1024x768 pixels is required to access SmartMonitoring.
- For your convenience, make a note of the serial numbers of your refrigerators in advance.
- A MyLiebherr customer profile is required to log in to the SmartMonitoring dashboard (registration possible during set-up).
- You will need additional accessories to set-up other appliances.
- Recommendation: one SmartCoolingHub per room
- If the distance between the SmartCoolingHub and the appliance is greater than 50 m, you should consult a qualified technician for installation.

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